

Resetting the X-ray Computer

-Aaron Sattler

11.4.2008

- First, make sure that goniometer is in its zeroed position
- Exit out of **APEX2**
- Exit out of **BIS**
- Exit out of all other programs
- Reset Computer
- When computer is booted up, put in password: *Bruker*
- First, open **BIS**
- Once open, confirm detector distance (*5.077cm*) and press okay
- Open **APEX2**
- In the toolbar menu, click on: *Instrument: Connect*
 - IP: 192.168.1.1
- Then in the toolbar menu, click on: *Instrument: Show Status*